# BENCH FUME HOOD

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HOTTES PARA BANCADA  
SORBONNES SUR LES PAILLASSES

## OVER BENCH

The range of Over Bench Fume Hoods is divided into three distinct models, the Cairo, the VWS and the PMMA Fume Hood, each of them has differentiating characteristics, sharing only the fact that they need a bench. Moreover, the materials, design and purpose differ that allows a more complete line, however these are simpler than the lines previously presented, not presenting, for example, the ability to adapt to the services that the user may need.

This range of Industrial Laborm fume hoods is developed according to standards:

**EN 14175-1: 2003** - Chemistry Fume Hoods - Part 1: Definition and Dimensions

**EN 14175-2: 2003** - Chemistry Fume Hoods - Part 2: Performance and Safety Requirements

**EN 14175-3: 2003** - Chemistry Fume Hoods - Part 3: Type Test Methods

**EN 13150: 2001** - Laboratory Benches - Dimensions, Safety Requirements and Test Methods.

**EN 61010-1: 2001** - Safety requirements for electrical equipment for measurement, control and use in the laboratory - Part 1: General requirements

**EN 61326-1: 2006** - Electrical equipment for measurement, control and use in the laboratory - Electromagnetic compatibility requirements - Part 1: General requirements.

CE marking according to the directives: **2014/30/EU** (Electromagnetic Compatibility) and **2014/35/EU** (low voltage).

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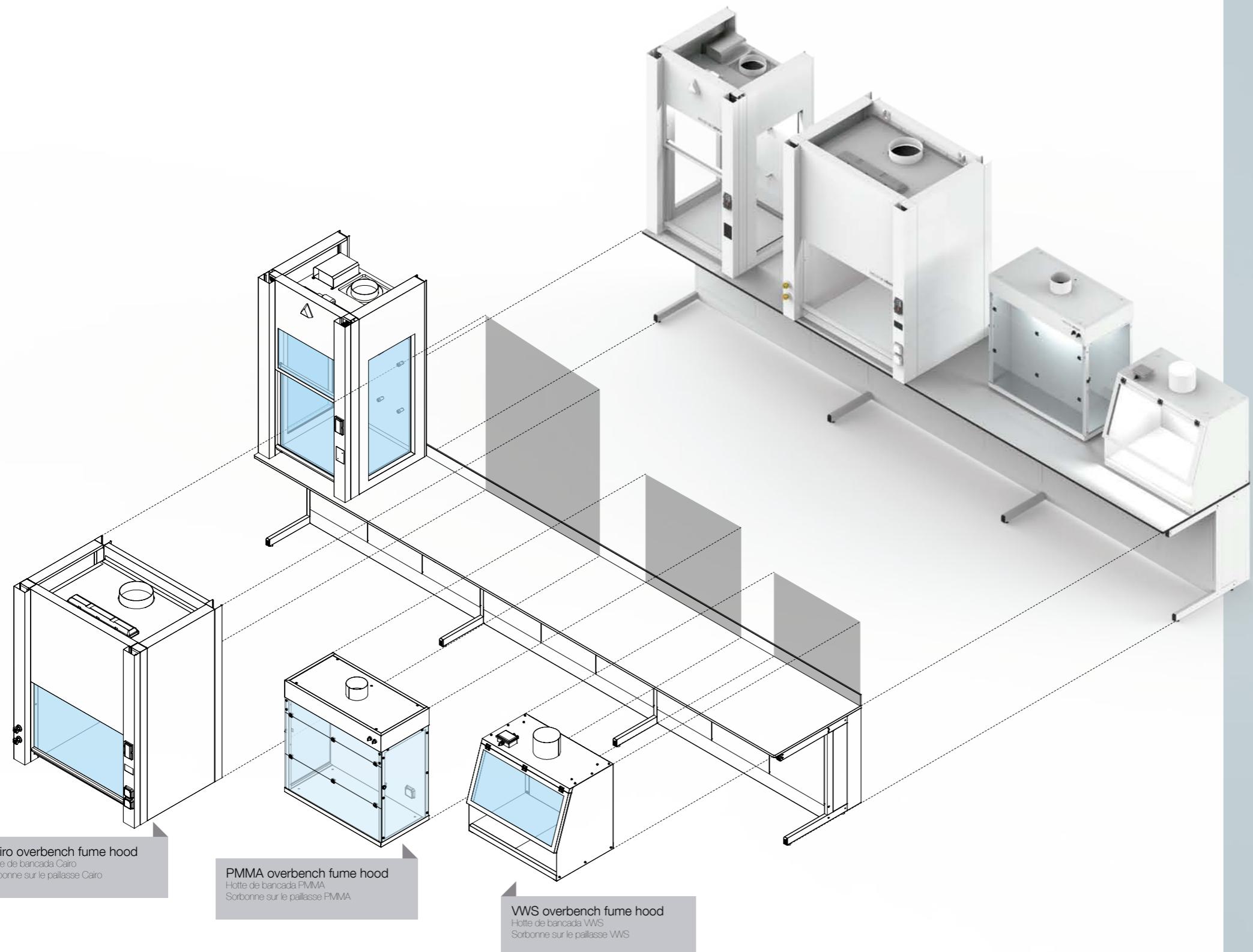
HOTTES CAIRO/VWS/PMMA  
SORBONNES CAIRO/VWS/PMMA

## CAIRO / VWS & PMMA

Cairo - Fume Hood designed for lab work that poses a high risk to users, is a fume hood consisting of two distinct parts, exterior body and inner lining. The exterior body serves as a support for the inner lining, is made of steel sheet powder coated with epoxy-polyester resins. It also supports the services that will be located on the sides. The inner lining is the part in direct contact with the chemical vapors, it is made of laminated compact. The standard model is supplied without any service, and is only supplied with a simple control CAV system, (Constant Air Volume).

VWS - Fume Hood developed for small laboratory work, made entirely of steel sheet powder coated with epoxy-polyester resins, this fume hood is divided into two parts, the body and a removable work surface. It is a Fume hood that does not allow a integration of services and is only provided with an ON/OFF switch that allows switching on and off the light and the extraction at the same time.

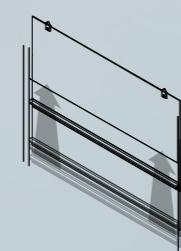
PMMA - Also referred to as an acrylic fume hood, it was designed for small laboratory work and it is divided into 4 parts. Structural body made of steel sheet powder coated with epoxy-polyester resins, serves as a support for the entire fume hood, it is in this part that we have the levelers that allow the work surface to be always level. Hinged front window with hinges with a maximum opening of 180°, can be manufactured in crystalline polymethyl methacrylate or in glass. Side walls, made of crystalline polymethyl methacrylate or steel sheet powder coated with epoxy-polyester resins. The fourth and last component of the fume hood is the work surface, and can be manufactured in stainless steel sheet or in solid ceramic stoneware. This hood allows the integration of electrical outlets in the inside and is provided with two ON/OFF switch that allows switching on and off the light and the extraction separately.

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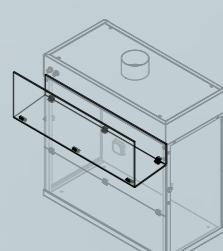
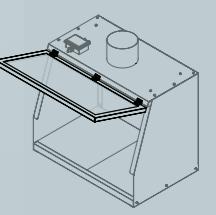
Fluxo de ar

Flux d'air

The Cairo Fume Hood comes with constant air volume (CAV), however can be equipped with VAV system (variable air volume), which improves the performance of the Fume Hood as well as its overall energy consumption.

In the same way the Fume Hood can also be equipped with automatic window closing system. The PMMA and VWS fume hoods only have an ON/OFF button for extraction and luminaire.

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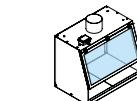
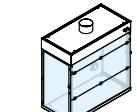
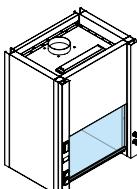
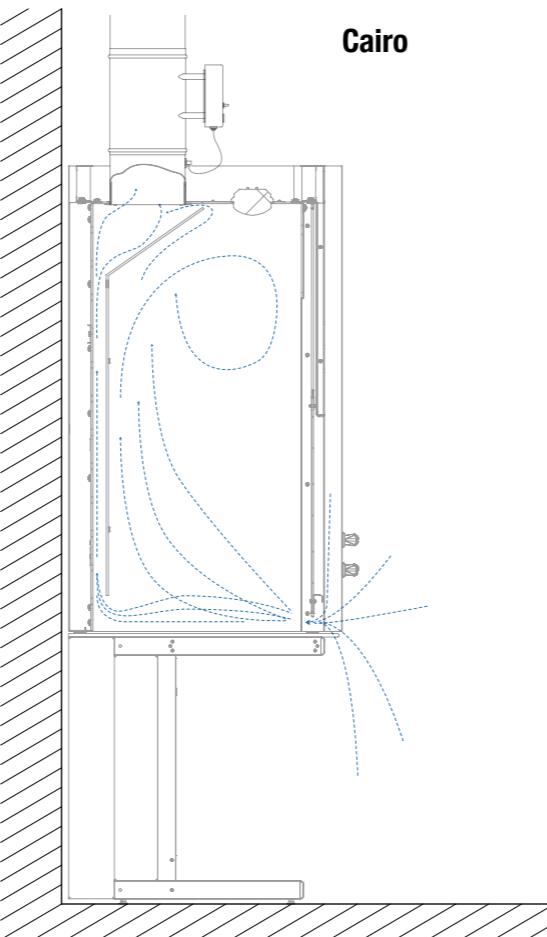
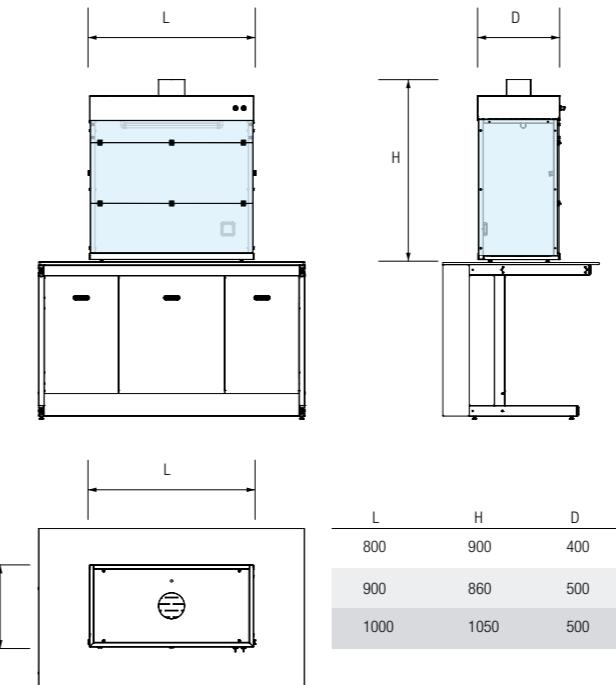
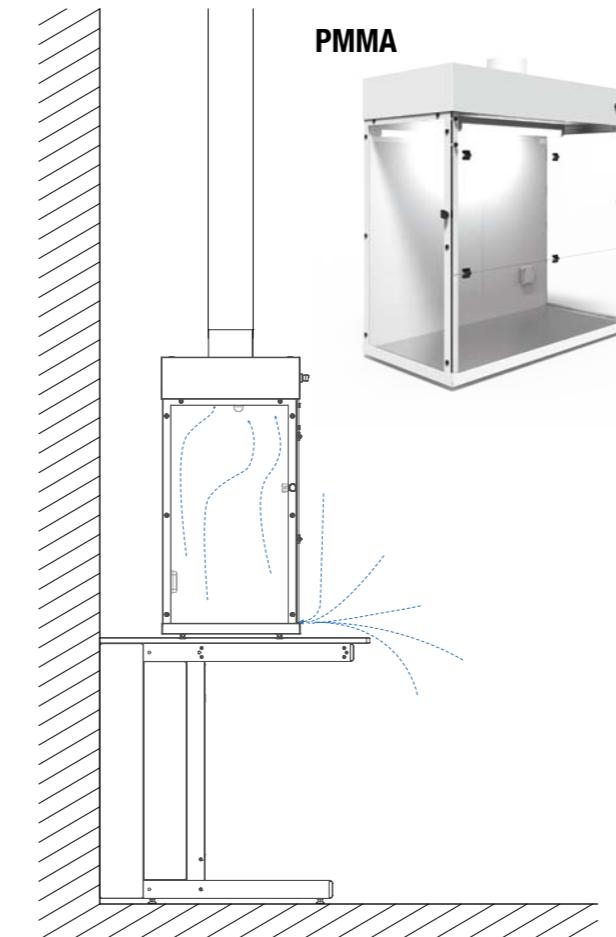
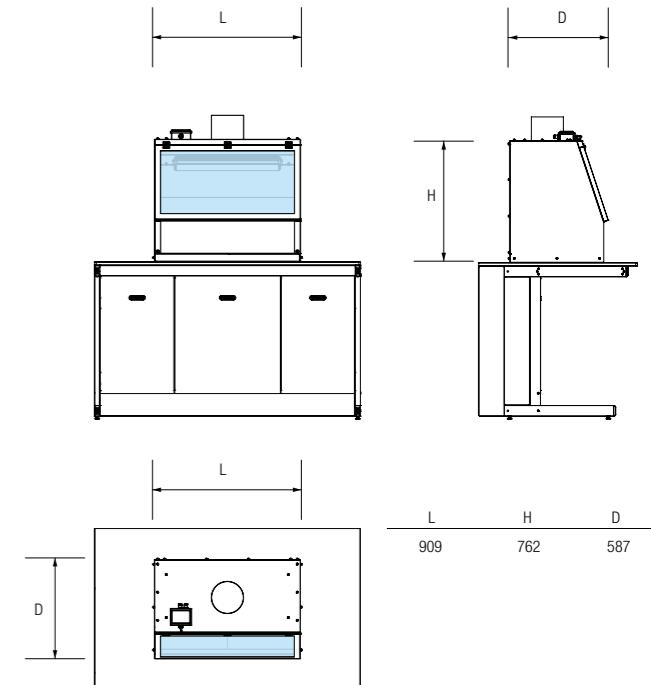
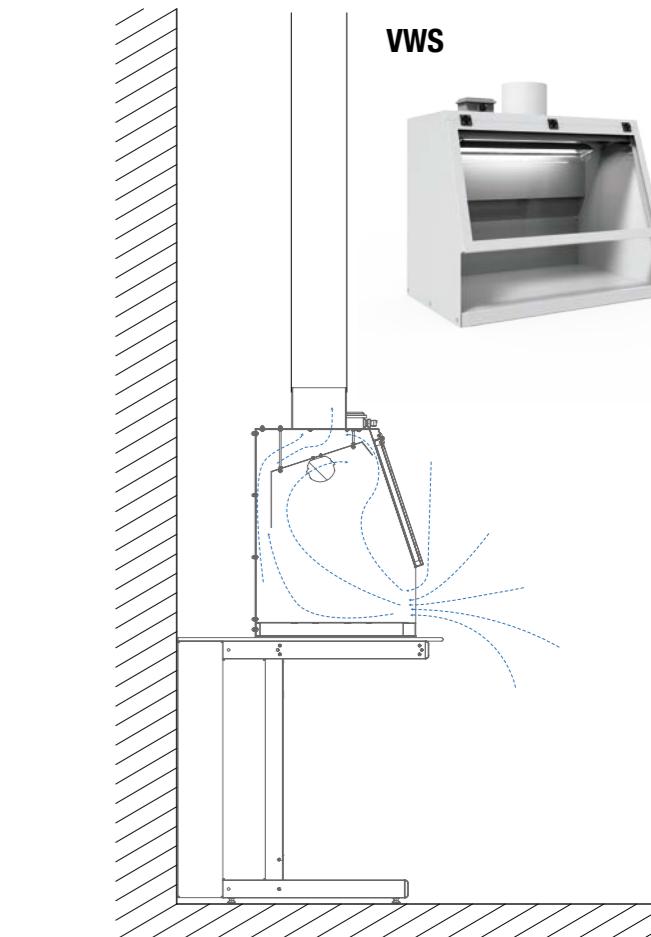
## Electrical information

Informação eléctrica  
Informations électriques

The ENS Fume Hood is a class I equipment, so all accessible metal parts are connected to the earth terminal, which must be connected to the earth protective circuit of the building. The terminal strips for electrically connecting the Fume Hood are located in the electrical cabinet on the control panels.

The equipment has all necessary magnetothermic protections, namely:

- Thermal protection of the fan (s);
- Protection circuit breaker of the circuit of the sockets;
- circuit breaker protection circuit breaker;
- Protection circuit breaker for the control circuit.

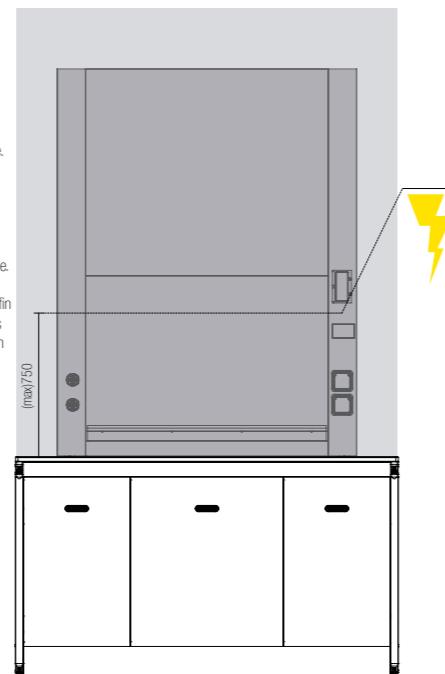
The installation site must have a circuit-breaker to ensure that in the case of other equipment connected to the circuit-breaker, these shall not be without power in case of maintenance of the Fume Hood.

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

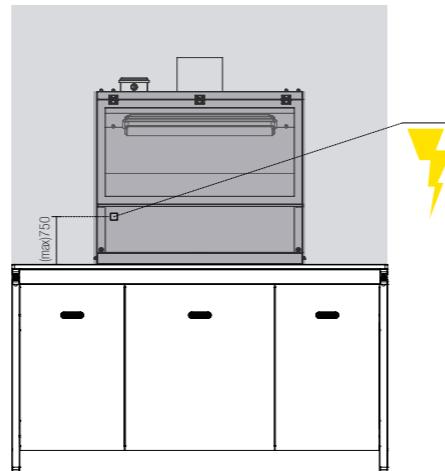
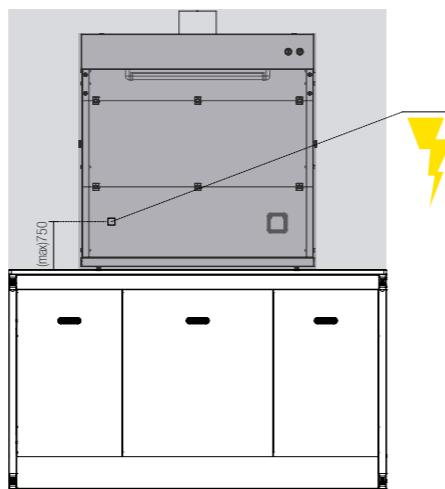
CAIRO

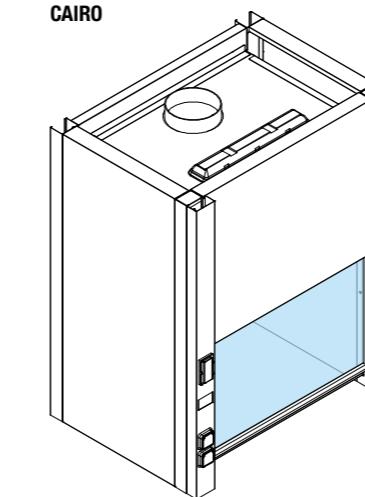
Cairo Fume Hood Standard for a widespread use of chemical material, with cold rolled sheet metal exterior and phenolic compact inner lining. This is the premium model of the line, presenting the possibility of incorporating VAV system and automatic sash closing. There is also a version with glazed sides and a central version with glazed sides, baffles and bottom in glass.

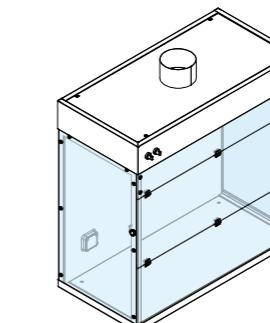PMMA

PMMA or Acrylic Fume Hood, designed for small jobs and small quantities of chemicals, with cold rolled steel sheet structure, stainless steel surface and sides and window in polymethyl methacrylate or glass. It's a very versatile fume hood, and is supplied with levelers which allows a work surface always level, does not present the possibilities of systems that the Cairo model presents, nevertheless it makes possible to have electrical services in the interior.

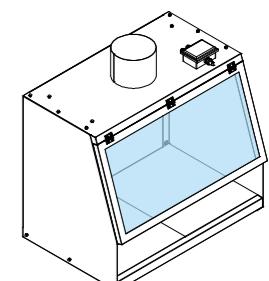VWS

VWS Fume Hood designed for small jobs and small quantities of chemicals. Built entirely of cold rolled steel sheet with a glazed window. It's a very versatile fume hood, however it does not present any possibility of integration of systems like Cairo or even of services like the PMMA, thus being the simplest fume hood that Industrial Laborum presents.

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