# **PATHOLOGY ANATOMY FUME HOOD**

HOTTE DE ANATOMIA PATOLÓGICA  
SORBONNE DE ANATOMIE PATHOLOGIQUE

# PATHOLOGIC ANATOMY

## FUME HOOD

HOTTE DE ANATOMIA PATOLÓGICA

SORBONNE D'ANATOMIE PATHOLOGIQUE

## PATHOLOGIC ANATOMY

The Pathologic Anatomy Fume hood is a safety equipment designed for laboratory work that poses a high risk to users. Their function is to control the user's exposure to gases, fumes, vapors, aerosols, dangerous particles and other work with volatile substances from the washing of materials, in chemistry laboratories and pathological anatomy laboratories.

This is the premium model of Industrial Laborum developed according to standards:

**EN 14175-1: 2003** - Chemistry Fume Hoods - Part 1: Definition and Dimensions

**EN 14175-2: 2003** - Chemistry Fume Hoods - Part 2: Performance and Safety Requirements

**EN 14175-3: 2003** - Chemistry Fume Hoods - Part 3: Type Test Methods

**EN 13150: 2001** - Laboratory Benches - Dimensions, Safety Requirements and Test Methods.

**EN 61010-1: 2001** - Safety requirements for electrical equipment for measurement, control and use in the laboratory - Part 1: General requirements

**EN 61326-1: 2006** - Electrical equipment for measurement, control and use in the laboratory - Electromagnetic compatibility requirements - Part 1: General requirements.

CE marking according to the directives: **2014/30/EU** (Electromagnetic Compatibility) and **2014/35/EU** (low voltage).

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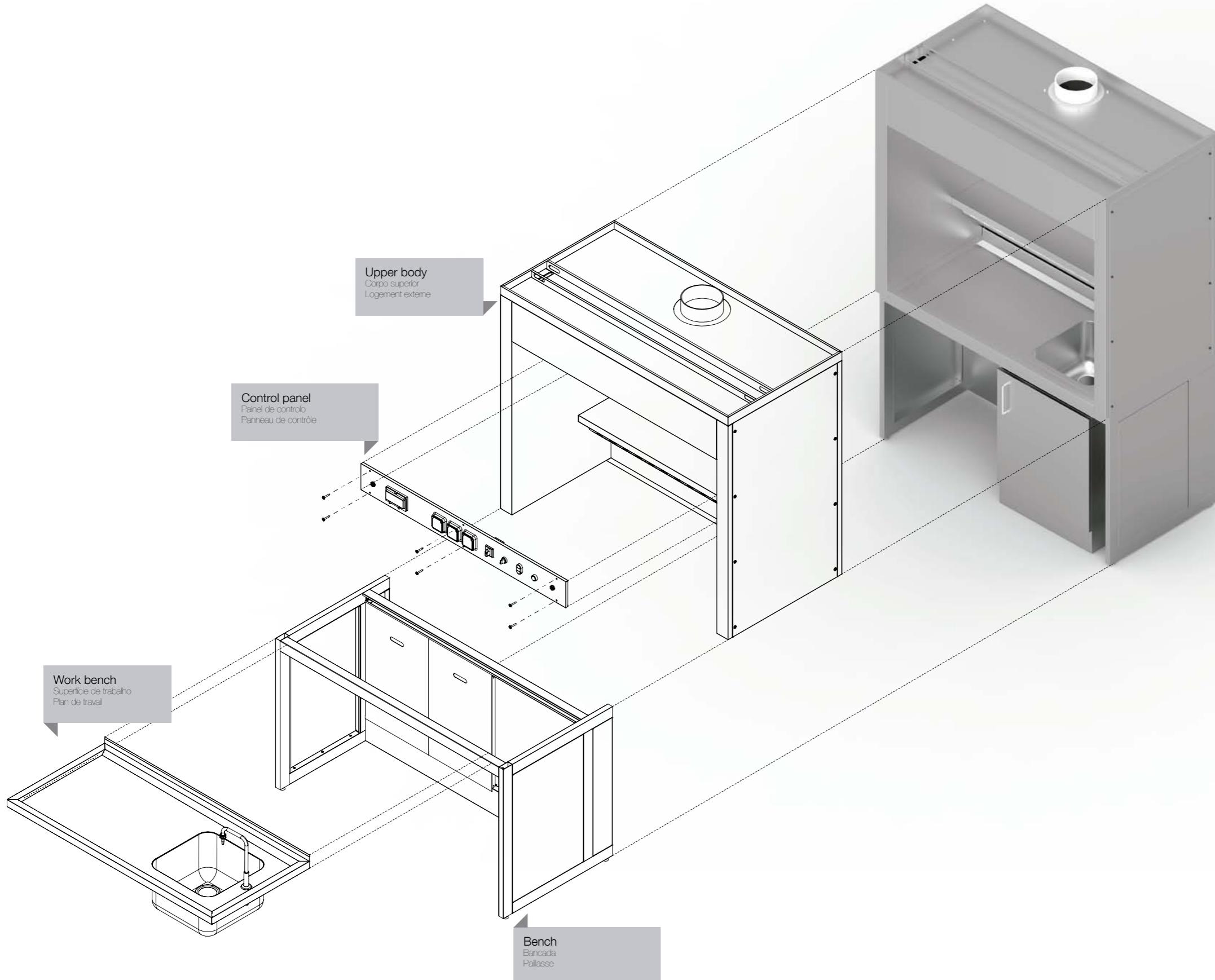
SORBONNE D'ANATOMIE PATHOLOGIQUE

## FUME HOOD

The Pathologic Anatomy Fume Hood is a confined space, accessible from the front. To maintain the containment inside the hood, a negative pressure is created in relation to the outside through the connection to the air extraction system.

The Pathologic Anatomy Fume Hood consists of:

- Workbench - The main function is to support the working surface and the upper body of the fume hood, built in tube and panels made of AISI 316 stainless steel sheet with high chemical resistance;
- Upper body - Fully built in AISI 316 austenitic stainless steel sheet, with high corrosion resistance. Connection of the various elements through screws, allowing easy disassembly and replacement of components. Interior of the upper body designed in order to favor the laminar flow of air, and to provide spaces for storage and placement of utensils.
- Working surface - Made of AISI 316 stainless steel sheet, it has a sink, a water faucet and a water jet system for horizontal washing of the working plane.

This fume hood has two extraction zones: a lower zone on the washing zone that allows suctioning at the work surface level, and an upper zone to remove vapors and more volatile gases.

Our Standard models are supplied with a simple CAV system (Constant Air Volume), and the services are a photoelectric cell water faucet, are also supplied with a sink on the worktop, as well as two or three IP55 single-phase sockets. The controls and electrical plugs are placed on the control panel above the first extraction zone, here also we find the photoelectric cells that drive the washing system of the work surface, as well as the crusher button, the control button of the flow extraction and the luminaire, finally has all the circuit breakers and protections of electrical components.

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
## Air flow

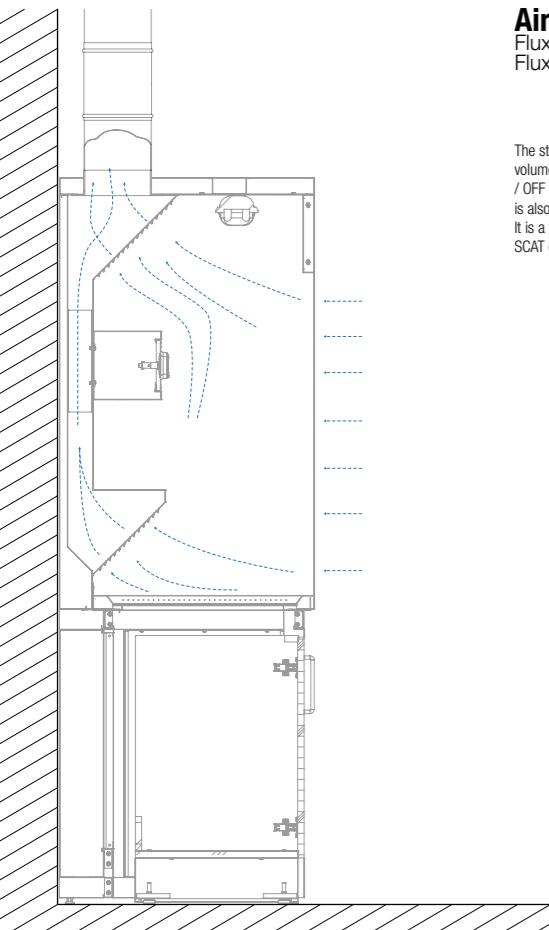Fluxo de ar  
Flux d'air

The standard Fume Hood comes with constant air volume (CAV), with the simplified system of an ON / OFF switch to turn the exhaust on and off, there is also a switch to turn the luminaire on and off. It is a fume hood that can be equipped with the SCAT chemical waste collection system.

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
## Standard services

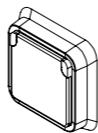Serviços standard  
Services standard

2/3 schuko-type outlets (type F),  
230V, protection class IP55\*\*

2/3 tomadas do tipo schuko (tipo F), 230V, classe de proteção IP55\*\*  
2/3 prises de type schuko (type F), 230V, degré de protection IP55 \*\*

One cold water valve with photoelectric cell activation

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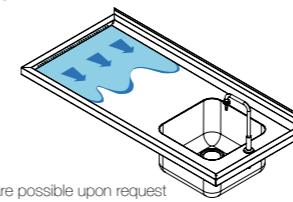
Work plan water jet cleaning system,  
activated by photoelectric cell

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\*\*model 1200 with two sockets and model 1500 and 1800 with 3 sockets, the location of the sockets will be on the front panels of the fume hood, and you must specify if you want in any other location.

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## Electrical information

Informação eléctrica  
Informations électriques

The Fume Hood is a class I equipment, so all accessible metal parts are connected to the earth terminal, which must be connected to the earth protective circuit of the building. The terminal strips for electrically connecting the Fume Hood are located in the electrical cabinet on the control panels. The equipment has all necessary magnetothermic protections, namely:

- Thermal protection of the fan(s);
- Protection circuit breaker of the circuit of the sockets;
- Circuit breaker protection of the lighting system;
- Protection circuit breaker for the control circuit.

The installation site must have a circuit-breaker to ensure that in the case of other equipment connected to the circuit-breaker, these shall not be without power in case of maintenance of the Fume Hood.

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## Control Panel

Painel de controlo

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Consisting of fan controls, electric protection cabinet, lighting and shredder pushbutton, photoelectric cells for tap and wash operation on the working plan.

The ventilation control has two indicators: Green indicator - Lit indicates that the fan is running.

Red flag - Lit indicates that the fan is off.

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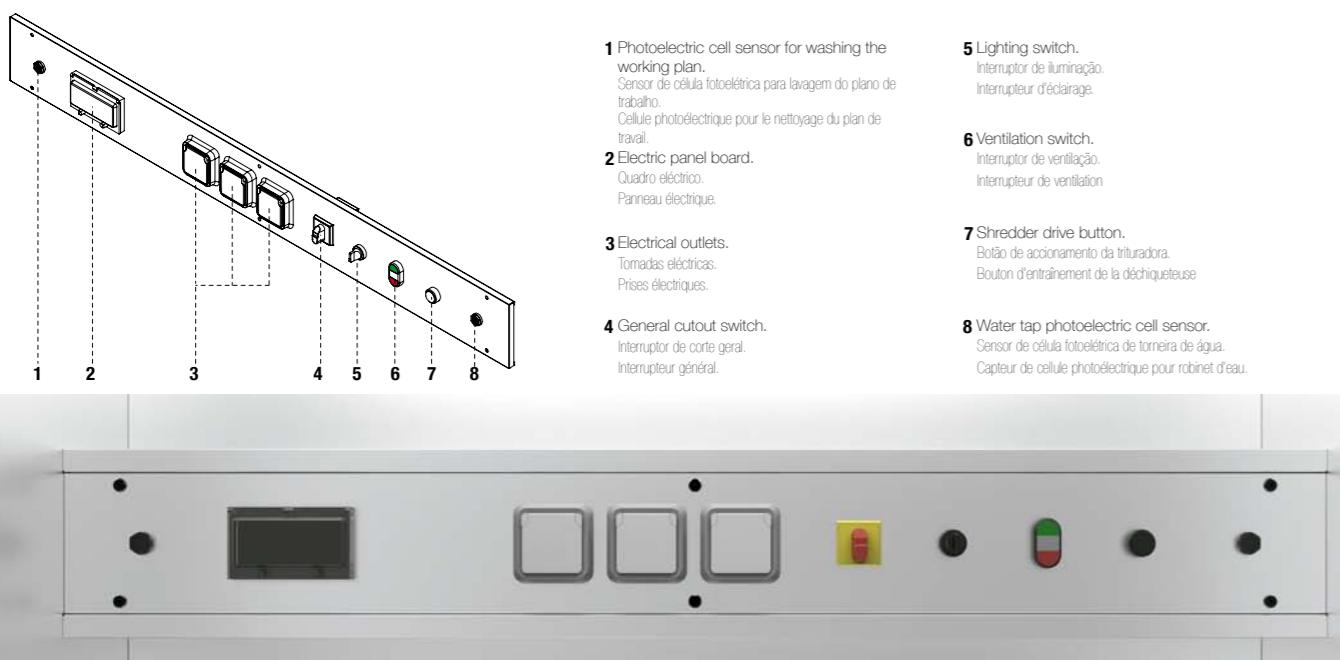
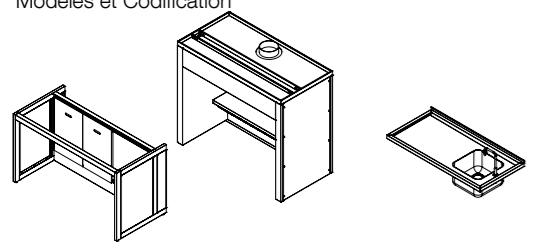
**STAINLESS STEEL**  
AÇO INOX  
ACIER INOXYDABLE

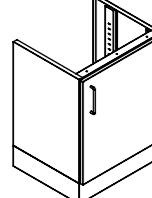
**STAINLESS STEEL**  
AÇO INOX  
ACIER INOXYDABLE

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