WASHING FUME HOOD

HOTTE DE LAVAGEM
SORBONNE DE LAVAGE

WASHING FUME HOOD
HOTTE DE LAVAGEM
SORBONNE DE LAVAGE

WASHING

The Washing Fume Hood is a safety equipment designed for laboratory work that poses a high risk to users. Their function is to control the user's exposure to gases, fumes, vapors, aerosols, dangerous particles and other work with volatile substances from the washing of materials, in chemistry laboratories and pathological anatomy laboratories.

This is the premium model of Industrial Laborum developed according to standards:
EN 14175-1: 2003 - Chemistry Fume Hood - Part 1: Definition and Dimensions
EN 14175-2: 2003 - Chemistry Fume Hood - Part 2: Performance and Safety Requirements
EN 14175-3: 2003 - Chemistry Fume Hood - Part 3: Type Test Methods
EN 13150: 2001 - Laboratory Benches - Dimensions, Safety Requirements and Test Methods.
EN 61010-1: 2001 - Safety requirements for electrical equipment for measurement, control and use in the laboratory - Part 1: General requirements
EN 61326-1: 2006 - Electrical equipment for measurement, control and use in the laboratory - Electromagnetic compatibility requirements - Part 1: General requirements.
CE marking according to the directives: **2014/30/EU** (Electromagnetic Compatibility) and **2014/35/EU** (low voltage).

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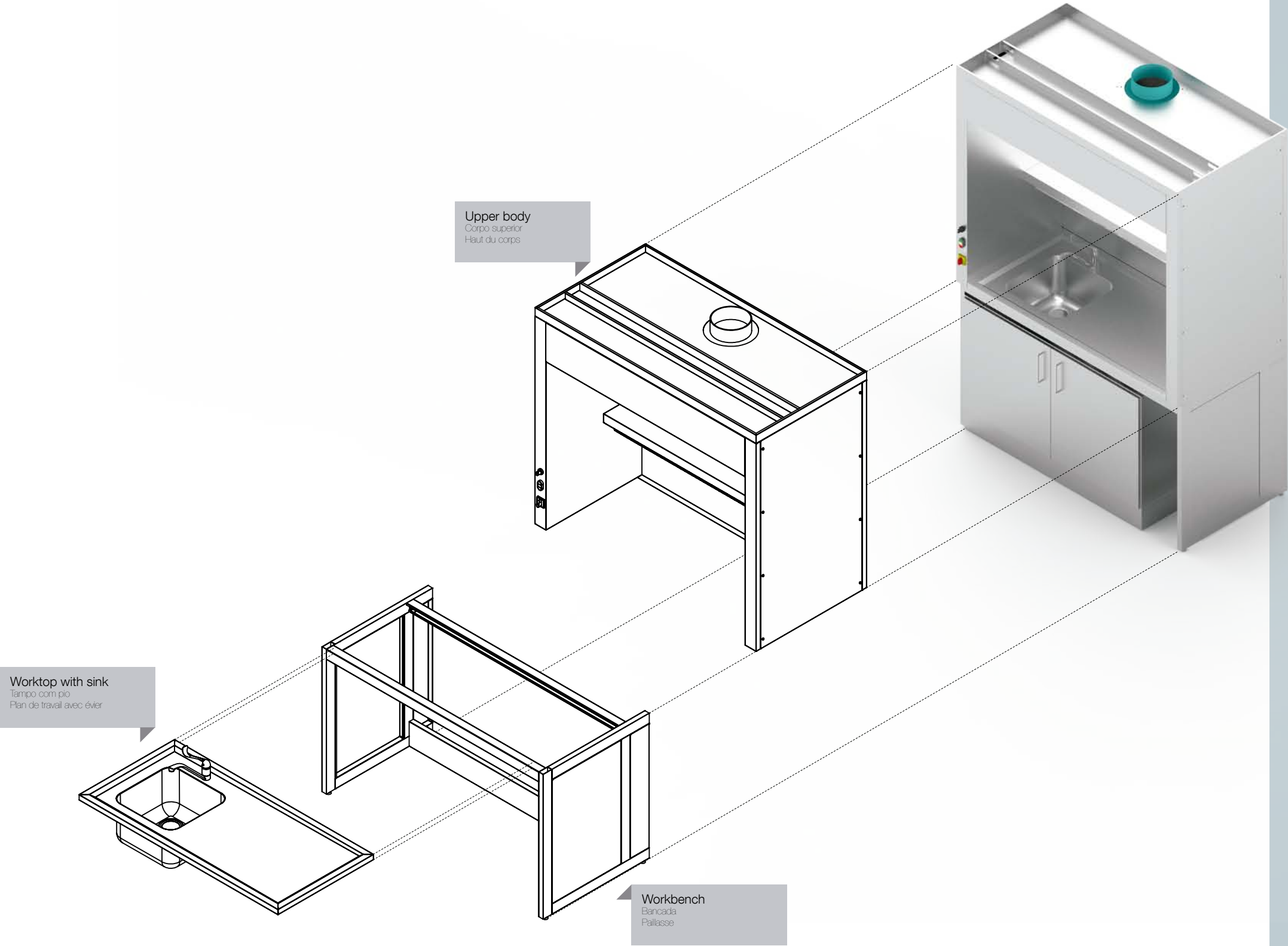
WASHING FUME HOOD
HOTTE DE LAVAGEM
SORBONNE DE LAVAGE

WASHING

The Washing Fume Hood work area is a confined space, accessible from the front. To maintain the containment inside the Fume Hood, a negative pressure is created in relation to the outside through the connection to the air extraction system.

- The washing fume hood consists of:
- Workbench - The main function of the workbench is the support to the work surface and to the upper body of the fume hood, built in tube and panels made of AISI 316 stainless steel sheet with high chemical resistance;
 - Upper body - Fully built in AISI 316 austenitic stainless steel plate with high corrosion resistance or with polypropylene interior. Connection of the various elements through screws, allowing easy disassembly and replacement of components. Interior of the upper body designed in order to favor the laminar flow of air, and to provide spaces for storage and placement of utensils;
 - Working surface - Made of AISI 316 stainless steel sheet, polypropylene or epoxy resin, it has a sink for washing and one or two water mixer taps.

This Fume Hood has two extraction zones: a lower zone on the washing zone that allows suctioning at the work surface level, and an upper zone to remove vapors and more volatile gases. Our Standard models come with a simple CAV (Constant Air Volume) system, and the services are a water mixing faucet plus one or two sinks on the top. The controls are placed on the left side of the fume hood, here is also the power cut button.

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Air flow
Fluxo de ar
Flux d'air

The standard Fume Hood comes with constant air volume (CAV), with the simplified system of an ON/OFF switch to turn the exhaust on and off, there is also a switch to turn the luminaire on and off. It is a fume hood that can be equipped with the SCAT chemical waste collection system.

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
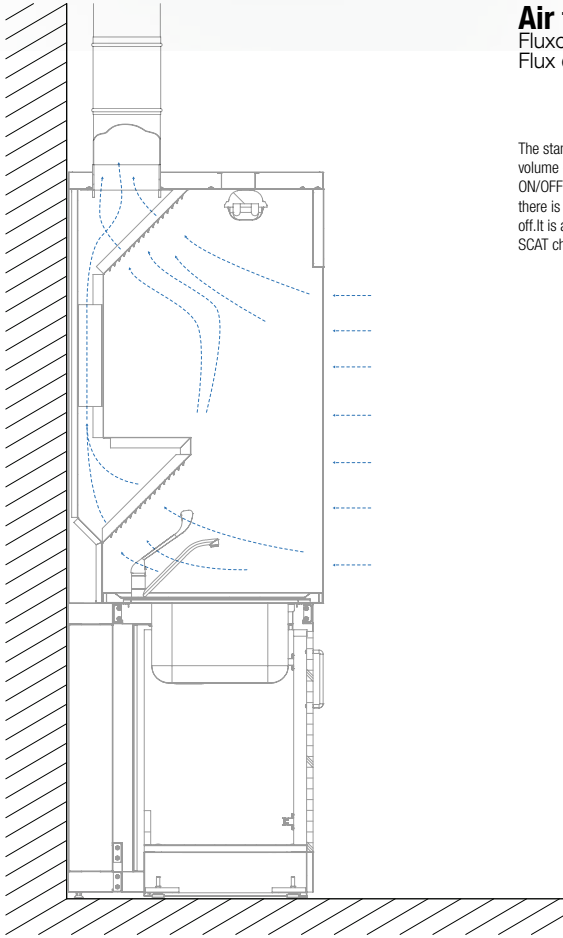
Table with 5 columns: Internal volume, Minimum flow rate, Nominal output diameter, Pressure drop. Rows for model sizes 1200, 1500, 1800.

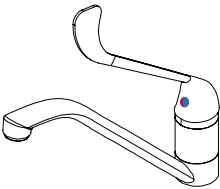Standard services
Serviços standard
Services standard

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Electrical information
Informação eléctrica
Informations électriques

The Fume Hood is a class I equipment, so all accessible metal parts are connected to the earth terminal, which must be connected to the earth protective circuit of the building. The terminal strips for electrically connecting the Fume Hood are located in the electrical cabinet on the control panels.
The equipment has all necessary magnetothermic protections, namely:
- Thermal protection of the fan(s);
- Disjuntor de proteção do circuito de iluminação;
- Circuit breaker protection for lighting system;
- Protection circuit breaker for the control circuit.


The installation site must have a circuit-breaker to ensure that in the case of other equipment connected to the circuit-breaker, these shall not be without power in case of maintenance of the Fume Hood.

Table with 5 columns: model size, Power Supply Voltage, Frequency, Power supply, Installed power. Rows for model sizes 1200, 1500, 1800.

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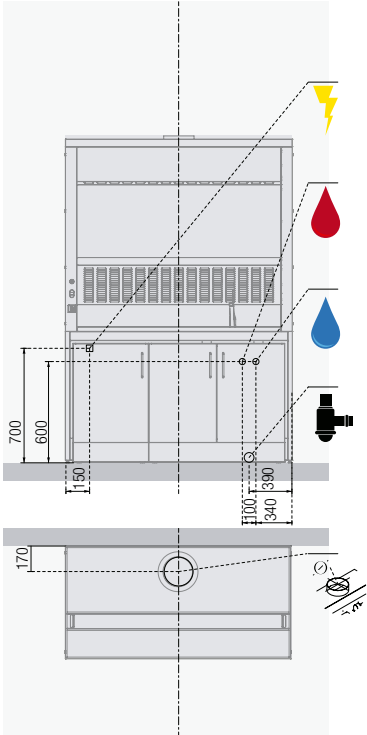
Services pre-installation
Pré-instalação de serviços
Pré-installation des services

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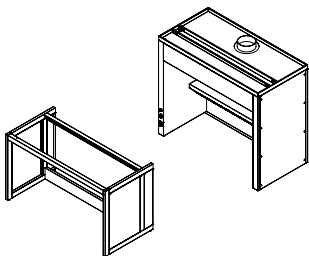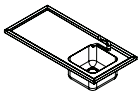
STAINLESS STEEL
AÇO INOX
ACIER INOXYDABLE


STAINLESS STEEL
AÇO INOX
ACIER INOXYDABLE

L x D x H*				
1200x750x2140	110-0001	110-1011	110-2001	
1500x750x2140	110-0002	110-1012	110-0003	
1800x750x2140	110-2004	110-1004	110-2005	110-0004

L x D x H*			
1200x750x2140	110-0021		
1500x750x2140	110-0022	110-0023	
1800x750x2140			110-0024

L x D x H*			
1200x750x2140	110-0011	110-2011	
1500x750x2140	110-0012	110-0013	
1800x750x2140	110-2014		110-0014

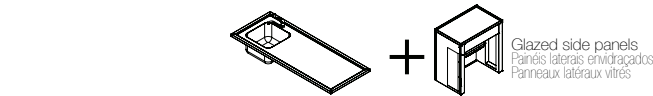
L x D x H*		
1200x750x2140	110-1001	
1500x750x2140	110-1002	
1800x750x2140		

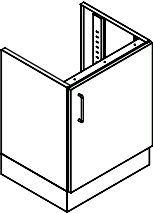*Length x Depth x Height
*Comprimento x Profundidade x Altura
*Longueur x Profondeur x Hauteur

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